

Rechargeable memory battery replacement


The internal battery has been found to have a shelf life of over 6 months. With normal use, the internal battery should not need to be replaced.

The symptom for a dead battery is no temperature display, when the Premier unit is turned off. If the temperature display comes on when the Premier is turned on, the battery may simply need re-charging, which takes about 8 hours of normal use. If after the re-charging time has elapsed and the temperature display is off when the Premier is turned off, please follow these instructions for its replacement or return the unit to Criminalistics, Inc. and we can replace the battery for you, for a nominal service charge.

- 1) Remove the 2 securing screws for the front panel.
- 2) Gently tilt the front panel forward, the AA NiMh battery will be visible with a securing tie wrap.



- 3) Cut the tie wrap, remove the AA battery and replace with either a NiCd or NiMH AA cell. Do not use an alkaline cell, except for emergency situations. Alkaline cells are not rechargeable.

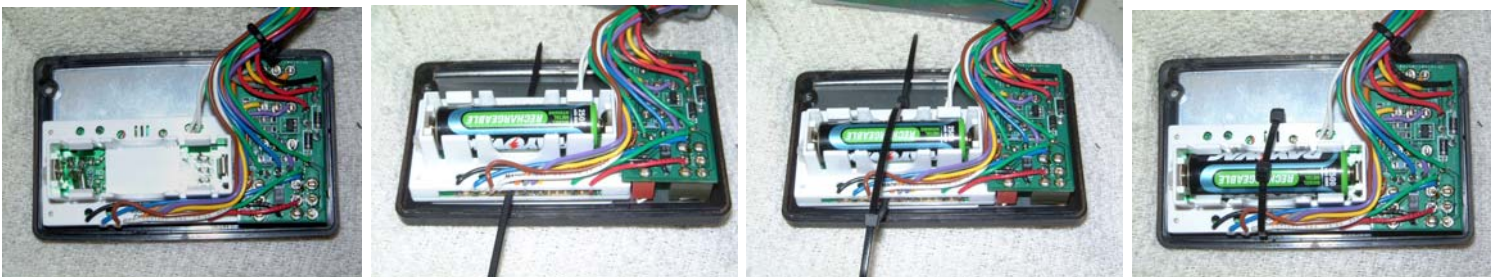

- 4) Replace the cut tie wrap (using 2 small tie wraps recommended) to secure the AA cell in place.
- 5) Carefully reseal and resecure the front panel to the unit.